

Identifier: Versa 20-MAN-0	Revision: 0	Effective Date: 10/7/15	 QUANTUM N · O · R · T · H · W · E · S · T
Web Page: http://www.qnw.com/versa-20-product-family/			
Description: Peltier-controlled cuvette holder with two optical ports for use in UV/Vis spectrophotometers			

QNW Versa 20 User's Guide

Quantum Northwest, Inc.
22910 E Appleway Avenue, Suite 4
Liberty Lake, WA 99019-8606
Phone: (509) 624-9290
Fax: (509) 624-9488
Service email: service@qnw.com
www.qnw.com

Temperature-Controlled Cuvette Holders
and Custom Instrumentation

Notices

Copyright

Copyright © 2015 Quantum Northwest, Inc. All rights reserved.

No part of this publication may be reproduced, transmitted, transcribed, stored in a retrieval system, or translated into any language, in any form or by any means, electronic, mechanical, photocopying, recording, or otherwise, without prior written permission from Quantum Northwest, Inc. All copyright, confidential information, patents, design rights and all other intellectual property rights of whatsoever nature contained herein are and shall remain the sole and exclusive property of Quantum Northwest, Inc. The information furnished herein is believed to be accurate and reliable. However, no responsibility is assumed by Quantum Northwest, Inc. for its use, or for any infringements of patents or other rights of third parties resulting from its use.

The Quantum Northwest name, QNW acronym and Quantum Northwest logo are trademarks or registered trademarks of Quantum Northwest, Inc. All other trademarks are the property of their respective owners.

Quantum Northwest products must only be used in the manner described in the Quantum Northwest produce user guides. Any other use may result in damage to the product or personal injury. Quantum Northwest is not responsible for any damages caused, in whole or in part, by improper use of the products, unauthorized alterations, adjustments or modifications to the products, failure to comply with procedures in Quantum Northwest product user guides, or use of the products in violation of applicable laws, rules or regulations. Quantum Northwest shall not be liable for errors or damages on products or performance due to using non-Quantum Northwest genuine parts.

Warranty

The material contained in this document is provided "as is," and is subject to being changed, without notice, in future editions. Further, to the

maximum extent permitted by applicable law, Quantum Northwest disclaims all warranties, either express or implied, with regard to this manual and any information contained herein, including but not limited to the implied warranties of merchantability and fitness for a particular purpose. Quantum Northwest shall not be liable for errors or for incidental or consequential damages in connection with the furnishing, use, or performance of this document or of any information contained herein. Should Quantum Northwest and the user have a separate written agreement with warranty terms covering the material in this document that conflict with these terms, the warranty terms in the separate agreement shall control.

Technology Licenses

The hardware and/or software described in this document are furnished under a license and may be used or copied only in accordance with the terms of such license.

Safety Notices

CAUTION

A CAUTION notice denotes a hazard. It calls attention to an operating procedure, practice, or the like that, if not correctly performed or adhered to, could result in damage to the product or loss of important data. Do not proceed beyond a CAUTION notice until the indicated conditions are fully understood and met.

WARNING

A WARNING notice denotes a hazard. It calls attention to an operating procedure, practice, or the like that, if not correctly performed or adhered to, could result in personal injury or death. Do not proceed beyond a WARNING notice until the indicated conditions are fully understood and met.

Table of Contents

1.	General Information	4
1.1.	User Documentation	4
1.2.	Modules, Covers and Panels	4
1.3.	Other Precautions	4
1.4.	Warning and Other Symbols	5
1.5.	CE Compliance	5
2.	Specifications	7
2.1.	Introduction: QNW Versa 20 Cuvette Holder with TC 1 Temperature Controller	7
2.2.	Measurement Category	8
2.3.	Pollution Degree	8
2.4.	Overvoltage Category	8
2.5.	Environmental Conditions	8
2.6.	Electrical Specifications	8
2.7.	Computer Requirements	9
3.	Installing, Starting and Stopping the QNW Versa 20	10
3.1.	Installing the QNW Versa 20	10
3.2.	Starting the QNW Versa 20	12
3.3.	Stopping the QNW Versa 20	13
4.	Step-by-Step Instructions for Using the QNW Versa 20 with TC 1 Temperature Controller	14
4.1.	Overview of TC 1 Temperature Controller	14
4.2.	Detailed Instructions	14
4.3.	External Computer Control	17
4.4.	Troubleshooting	17
4.5.	Error Messages	17
5.	Maintenance and Spare Parts	19
5.1.	Maintenance	19
5.2.	Spare Parts	19
	Appendix – Serial Communications for the TC 1 Temperature Controller	20

1. General Information

Thank you for purchasing a Quantum Northwest **Versa 20**. We want you to enjoy many years of faithful service from your instrument. If you have any questions, feel free to contact us directly through our website, www.qnw.com/contact-us/, or by email at service@qnw.com.

Your Quantum Northwest cuvette holder and accessories have been carefully designed so that when used properly, you have an accurate, fast, flexible and safe temperature control system.

Information about safety practices appears throughout the documentation provided with your instrument and accessories to help you safely operate the instrument and accessories. Before using the instrument or accessories, you must thoroughly read these safety practices. ALWAYS operate the instrument and accessories in accordance with these safety practices.

1.1. User Documentation

You have been provided with the following documentation to help you set up and operate your Quantum Northwest **Versa 20** cuvette holder:

- This manual, with safety practices and hazards information, instructions for installing and maintaining the components of the QNW **Versa 20**, and troubleshooting information.
- The **T-App** Help file, located on the **T-App** CD, if purchased (Quantum Northwest>**T-App** Help).

1.2. Modules, Covers and Panels

The spectrophotometer module that operators and other personnel will routinely access is the instrument's sample compartment, which is designed to contain the QNW **Versa 20**. Customers may insert cuvettes and cuvette accessories into the QNW **Versa 20** while the sample compartment is open.

Any other panels or covers that are retained by screws on the cuvette holder and accessories may be opened ONLY by Quantum Northwest service technicians.

1.3. Other Precautions

Do not block any ventilation grills present on the QNW **TC 1** Temperature Controller.

Use of the QNW **Versa 20** system and accessories may involve materials, solvents and solutions that are flammable, corrosive, toxic or otherwise hazardous. Careless, improper, or unskilled use of such materials, solvents and solutions can create explosion hazards, fire hazards, toxicity and other hazards which can result in death, serious personal injury, and damage to equipment and property.

ALWAYS ensure that laboratory safety practices governing the use, handling and disposal of such materials are strictly observed. These safety practices should include the wearing of appropriate safety clothing and safety glasses.

1.4. Warning and Other Symbols

The following is a list of symbols that appear in conjunction with warnings on the QNW cuvette holder and accessories. The hazard they describe is also shown.

A triangular symbol indicates a warning. The meanings of the symbols that may appear alongside warnings in the documentation or on the instrument itself are as follows.

	<p>When used on a warning label, the user is in danger of an electrical shock.</p>
	<p>When used on warning labels attached to the instrument, refer to the relevant operation or service manual for the correct procedure referred to by that warning label.</p>
	<p>When attached to the rear of the instrument, indicates that the product complies with the requirements of one or more EU directives.</p>

1.5. CE Compliance

Your QNW **Versa 20** cuvette holder has been designed to comply with the requirements of the Electromagnetic Compatibility (EMC) Directive and the Low Voltage (electrical safety) Directive (commonly referred to as the LVD) of the European Union. Quantum Northwest has confirmed that each product complies with the relevant Directives by testing a prototype against the prescribed EN (European Norm) standards.

Proof that a product complies with these directives is indicated by:

- the CE Marking appearing on the rear of the product, and
- the following Declaration of Conformity. The Declaration of Conformity is the legal declaration by Quantum Northwest that the product complies with the directives listed above, and shows the EN standards to which the product was tested to demonstrate compliance.

EC Declaration of Conformity

We: Quantum Northwest, Inc.
Of: 22910 E Appleway Ave, Suite 4
Liberty Lake, WA 99019-8606
USA

Declare under our sole responsibility that the products:

Models: **TC 1**, **turret 6**, **t2x2 Sport**, **BATH 10**, External Temperature Probe

Types: Temperature Controller, Temperature-Controlled Cuvette Holder, Submersible Pump, Temperature Probe

Are in conformity with the following standards and relevant EC directives:

Emissions: EN 61326-1 (2013)

Immunity: EN 61326-1 (2013)

EU Directive: 2004/108/EC, EMC

I hereby declare that the equipment models named above have been assessed and found to comply with the relevant sections of the above-referenced specifications. The units comply with the relevant requirements of the applicable Legislation, and I am the person authorized to compile the technical documentation.

Signed by: 
Name: Enoch W. Small, Ph.D.
Position: President, Quantum Northwest, Inc.
Date: August 22, 2014
Location: Liberty Lake, Washington, USA

2. Specifications

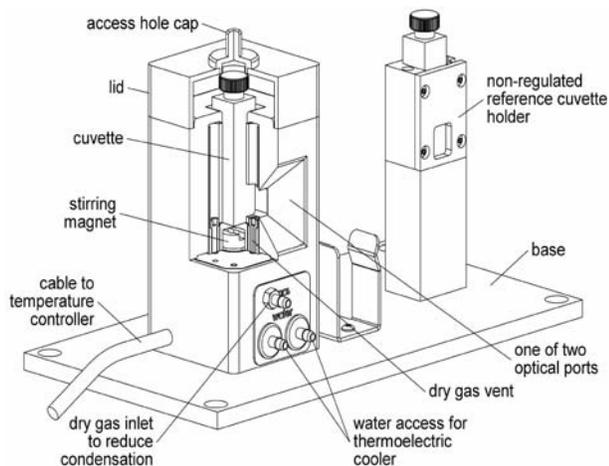


Figure 1. A typical configuration of the **Versa 20** for use in the Cary 300.



Figure 2. TC 1 Temperature Controller

2.1. Introduction: QNW Versa 20 Cuvette Holder with TC 1 Temperature Controller

Intended Use

The QNW **Versa 20** Cuvette Holder is designed to be inserted into a UV/Vis **spectrophotometer**, replacing the sample cuvette holder originally supplied with the instrument. The **Versa 20** has two optical ports for absorbance measurements. Temperature control is achieved using the **TC 1** Temperature Controller alone or along with a computer interfaced to the **TC 1** Temperature Controller box, using the QNW **T-App** software provided.

Specifications

- | | |
|-------------------------------------|------------------------------|
| • Temperature range | -15 to +110 °C* |
| • Temperature precision | ±0.01 °C |
| • Temperature accuracy | ± 0.2 °C from 0 °C to +80 °C |
| • Temperature reproducibility | better than ±0.05 °C |
| • Cuvette size (outside dimensions) | 12.5 x 12.5 mm |
| • Optical port dimensions | 12 mm high x 10 mm wide |
| • Probes accepted | Series 400 or Series 500 |
| • Cuvette z-height | 8.5 or 15 mm |

*Work below -15°C requires added windows and pre-cooled circulating fluid.

Note: The “z-height” of a cuvette is the distance between the bottom surface of the cuvette and the designed position for the optical center line, where the incident beam of light strikes the cuvette.

2.2. Measurement Category

The measurement category is IEC 61010:I. Do not to use this equipment for measurements within measurement categories II, III and IV.

2.3. Pollution Degree

The pollution degree is IEC61010:2. Pollution degree “2” applies to a normal indoor atmosphere.

2.4. Overvoltage Category

The overvoltage category (installation category) is CAT II. See IEC 664 & IEC 61010.

2.5. Environmental Conditions

The area should have a dust-free atmosphere with minimal drafts, vibrations, and corrosive fumes. For optimum performance, the ambient air is recommended to be between 20 and 25 °C, but can be from 5 to 40 °C. Relative humidity should be less than 80%. The instrument is designed for operation at 2,000 m elevation or less.

To avoid damage through spillage of solutions and samples being analyzed, the worktops should be covered with a material that is corrosion resistant and impervious to liquids.

Allow at least two inches of space on both sides, and six inches at the rear of the instruments to permit free air circulation. The power cord and all other connections are located at the rear of the instruments. The power switch is located on the front panel.

2.6. Electrical Specifications

Mains Supply

A standard power cable is provided based on the user's country requirements (18 AWG/115 V AC US/Canada; 1.0 mm/220 V AC international). The required supply voltage is 100-240 V AC (frequency 50 or 60 Hz). The mains voltage fluctuations must not exceed 10% of the nominal voltage.

The installation of electrical power supplies must comply with the rules and/or regulations imposed by local authorities responsible for the supply of electrical energy to the workplace.

All power supplies for the **TC 1** Temperature Controller must be single-phase, AC voltage, three-wire system (active, neutral, earth) and should be terminated at an appropriate power outlet receptacle that is within reach of the power cord. For safety reasons, a separate power outlet receptacle should be provided for each unit in the system. The use of extension cords or outlet adaptors is not recommended.

WARNING

Shock Hazard - Danger of electrocution. Good electrical grounding is essential to avoid potentially serious shock hazards. A 3-wire outlet with ground connection must be provided for the **Versa 20**. Make certain that power outlets are earth-grounded at the grounding pin.

CAUTION

Caution - Never connect or disconnect any cables while the **Versa 20/TC 1** Temperature Controller is switched on. Damage to the printed circuit boards may occur.

2.7. Computer Requirements

The **T-App** temperature application software used for external control of Quantum Northwest temperature controllers requires a personal computer using a Microsoft Windows™ operating system (XP, Vista, 7, or 8). The **T-App** software is normally provided on a CD, requiring a CD player be part of the computer. However, **T-App** may also be downloaded *via* the Internet from a QNW ftp site. Contact QNW should you wish to obtain the application by Internet.

3. Installing, Starting and Stopping the QNW Versa 20

3.1. Installing the QNW Versa 20

Unpacking

After accepting delivery, take the equipment to the installation site. Quantum Northwest instruments are inherently robust, and the packaging is designed to prevent internal damage. However, the contents form part of a precision measuring system and all packages should be handled with care. In transit, sharp jolts must be avoided and the packages should not be inverted or tilted unnecessarily. Markings on the shipping cartons generally indicate which side of the package should be kept on top.

Note: The **Versa 20** may have been already installed in the spectrophotometer by the spectrophotometer manufacturer.

Unpacking the equipment is your responsibility. As the packages are opened, ensure that you have received everything you ordered. If there are any discrepancies, notify the supplier. If any items are found to be damaged, immediately notify the carrier and supplier.

You should have received:

- 1x **Versa 20** Cuvette Holder with base plate
- 1x opaque lid
- 1x **TC 1** Temperature Controller
- 1x submersible pump
- 1x bucket for submersible pump
- 1x power cable
- 1x USB cable
- 1x CD containing **T-App** software (if ordered)
- 1x magnetic stir bar
- 1x vinyl tubing to connect water and gas
- This manual

Note: Any additional accessories ordered are not listed here.

Any differences from the original order should be referred immediately to your Quantum Northwest sales office. Do not discard any packaging components or filler materials.

Installing the QNW T-App Software (if ordered)

1. Insert the **T-App** program CD into the computer's CD-ROM drive. If the installation does not start automatically, locate the Setup.exe file in the root directory and run it. The installation process starts with a small black window that is shown during installation of the drivers needed to control the **Versa 20** through a USB connection. This window will then be replaced by the software installation window. Follow the onscreen prompts to complete the installation.
2. Connect the **Versa 20** to your computer using the USB cable provided. The New Hardware installation process will begin automatically and take a few moments to finish.

Installing the QNW Versa 20 Hardware



Figure 3. Submersible pump, bucket, electrical cable, and vinyl tubing for the **Versa 20**

1. Remove the existing sample compartment from the spectrophotometer. Install the **Versa 20** into the sample compartment and secure it in place.

Note: The **Versa 20** is supplied completely assembled. It is necessary only to make the appropriate connections.

2. Position the bucket to within 75 cm (30 in) above or below the spectrophotometer.

CAUTION



Caution – Position the bucket in a location that minimizes the risk of spilling the liquid contents.

3. **OPTIONAL** - If a gas such as nitrogen is to be used as dry gas for minimizing condensation at low temperatures, secure the tank to a solid object in the vicinity of the spectrophotometer.

CAUTION

Caution – When using dry gas do prevent condensation on the cuvette, ensure that the working area is adequately ventilated.

4. Cut the vinyl tubing provided into two lengths (three if you will be using the dry gas purge). Attach one length of tubing from the hose barb of the submersible pump, shown in Fig. 3, to either one of the water hose barbs on the **Versa 20**, labeled “water access for thermoelectric cooler” in Fig. 1. Attach another piece of tubing from the second **Versa 20** water hose barb to the bucket to serve as a drain tube. Add water to the bucket and lower the pump into the bucket, making sure that the pump is below the water line. Using the electrical cable from the pump, connect the pump to the socket on the back of the **TC 1** Temperature Controller. The pump will start automatically when the **TC 1** power is turned on later.
5. OPTIONAL - If you plan to work at low temperatures, connect a source of dry gas (typically nitrogen) to release upon the cuvette windows. Attach a length of tubing with $\frac{1}{8}$ -inch inside diameter to the small dry gas inlet hose barb on the base of the **Versa 20** (see Figure 1). Set the dry gas flow rate to about 50 cc/min.
6. Connect the **Versa 20** electrical cable to the 15-pin, VGA-type connector labeled “Sample” on the back of the **TC 1** Temperature Controller.
7. Plug the power cable into the back of the **TC 1** Temperature Controller and into a wall socket.

3.2. Starting the QNW Versa 20

1. Place the magnetic stir bar in a standard 1 x 1-cm square cuvette. Insert a liquid sample into the cuvette.
2. Place the cuvette in the **Versa 20**.
3. Place the plastic lid on the cuvette holder.
4. OPTIONAL - If you wish to monitor the temperature inside the cuvette, plug a standard Series 400 or Series 500 thermistor probe (not provided) into the $\frac{1}{4}$ - inch phone jack labeled “probe” in the back panel of the **TC 1**. Place the end of the probe in a region of the sample in the cuvette where it will not occlude the spectrometer light beam.
5. Turn on the **TC 1** controller using the switch on the back panel.¹

¹ When temperature control is turned on (see section 4.2), the **TC 1** controller supplies 12 V to the submersible pump, which will begin circulating water through the **Versa 20**. At this time, check for water leaks.

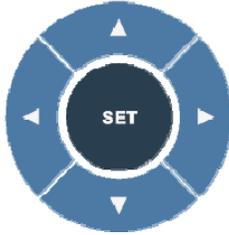


Figure 4. TC 1 Menu Button

6. **Brief Instructions:** The controller display begins on the **Display** page. Use the left or right arrows on the Menu Button (Figure 4) to cycle through the four pages of options, **Display** (the default screen), **Set Target Temperature**, **Set Stirring**, and **Set Ramping**. For example, to set a target temperature, press the right arrow once to go to the **Set Target Temperature** page. Once there, use the up and down arrows to set a target temperature, e.g. 37.0 °C. Press **SET** to initiate temperature control. While the temperature is changing, the green LED light above the display will flash slowly; when the temperature is stable, the LED will be steadily illuminated. To control stirring, go to the **Set Stirring** page. Use the up and down arrows to set a speed, e.g. 1200 rpm, and press **SET** to start the stirrer. **More details on the Menu Button are provided in Section 4.2 below.**

3.3. Stopping the QNW Versa 20

1. On the **Set Target Temperature** and **Set Stirring** pages, pressing the **SET** button without first making changes in numerical parameters will toggle the function ON/OFF. Check the **Display** page afterwards to confirm the system's status.
2. To completely turn off the **Versa 20**, turn off the power on the back of the **TC 1** controller. If a separate coolant pump was used, turn it off.

4. Step-by-Step Instructions for Using the QNW Versa 20 with TC 1 Temperature Controller

4.1. Overview of TC 1 Temperature Controller

The **TC 1 Temperature Controller**, shown in Figure 2, may be used in a stand-alone manner together with the **Versa 20** Peltier-controlled cuvette holder, or under PC control using QNW's **T-App** software or other vendors' software.

T-App is a program that permits external computer control of the **TC 1 Temperature Controller**. The temperature of the sample holder and a temperature sensed by an external probe may be plotted vs time. Simple text scripts may be used to automate multiple operations. **T-App** is provided with a USB cable to connect your computer to the **TC 1** drivers. Full instructions for using **T-App** may be found in the associated Help file.

4.2. Detailed Instructions

Using the Menu Button control arrows

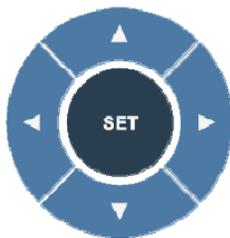


Figure 5. TC 1 Menu Button control arrows

Use the right and left arrows to cycle through four pages:

– Display – Set Target Temperature – Set Stirring – Set Ramping

Display

Holder = 37.0 °C Target = 37.0 °C Probe = 36.9 °C Ramp On Stir On

Figure 6. Example **TC 1** Display page

Display: This main page shows the actual cuvette *Holder* temperature, the *Target* temperature and a *Probe* temperature (if a probe is present). The page also shows whether the *Ramp* and *Stir* functions are on or off. After a few seconds of inactivity, all other pages will revert back to the **Display** page.

When seeking a temperature, the green LED on the front panel will flash slowly. When it has locked onto a new target temperature, the green LED will remain lit. A rapidly flashing red LED usually means a loose electrical cable or inadequate water flow for the Peltier unit.

Note: Pressing the **SET** button while the **Display** page is active results in no action. To start or stop controller functions, access the four other pages, **Set Target Temperature – Set Stirring – Set Ramping – Set Position**.

Set Target Temperature


Set Target Temp. Target = 37.0 °C Current= 23.6 °C Off
--

Figure 7. **TC 1** Set Temperature page

Set Target Temperature: To set the *Target* temperature, use the up and down arrows. Press **SET** to retain this new *Target* and initiate temperature control. The green LED light will begin flashing slowly as the device seeks the targeted temperature, and the page window will show the *Current* to be *ON*.

Discontinue Temperature Control: Press **SET** while the green LED light is lit or slowly flashing to discontinue temperature control. The LED light will turn off completely, and the page window will show the *Current* to be *Off*.

Set Stirring

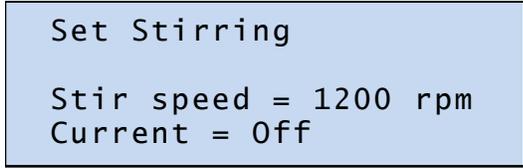


```
Set Stirring
Stir speed = 1200 rpm
Current = Off
```

Figure 8. TC 1 Set Stirring page

Set Stirring: To turn on magnetic *Stirring*, use the up and down arrows to choose an approximate stirring speed between 1 and 2500 rpm. Press **SET** to set the speed and initiate stirring. The page will update to show the *Current* value to be the same as *Stir speed*.

Discontinue Stirring: Press **SET** while the *Stirring* is on (shown by nonzero *Current* rpm) to turn the stirrer off. The page will update to show the *Current* speed to be *Off*.

Set Ramping

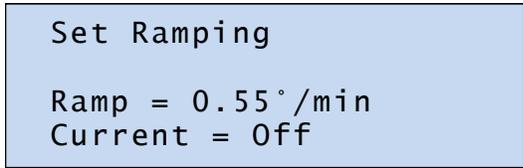

```
Set Ramping
Ramp = 0.55 °/min
Current = Off
```

Figure 9. TC 1 Set Ramping page

Set Ramping: To perform a temperature ramp, set the *Ramp* rate using the up and down arrows, and press **SET**. The page will update to show the *Current* rate in °/min to be the same as the *Ramp* rate. With ramping set, turning on temperature control will generate a linear ramp to the target temperature. When the sample holder reaches the target temperature, no further temperature change occurs, although the *Current* rate will remain as set.

The fastest possible ramp is determined by how fast the cuvette holder could reach the target temperature without ramping. Attempting to ramp too quickly, especially at high and low temperature extremes, will result in a nonlinear ramp. The slowest ramp that may be set on this page is 0.01 °C/minute. (If needed, much slower ramps may be set through software commands.)

Discontinue Ramping: Press **SET** while the *Ramp* is on (shown by nonzero *Current* °/min) to turn the *Ramp* off. The page will update to show the *Current* rate to be *Off*.

4.3. External Computer Control

All functions may be accessed either through Serial (RS 232) control or a USB located on the back of the **TC 1** Temperature Controller. You may write your own program or purchase the QNW application **T-App**. **T-App** will plot temperatures of the probe, cuvette tower, and even the Peltier element heat exchanger vs time. It will also permit you to set up temperature ramps. If you wish to do your own programming, please see the Appendix for communication instructions and the set of text commands that may be used and responses to the commands.

4.4. Troubleshooting

If the submersible pump has operated without being immersed in water, it may not prime properly. If this occurs, place one drop of detergent on the pump intake and try again.

If the **TC 1** display indicates that there is inadequate coolant, it may be that the circulating water is too warm. Add ice to the water to cool it down.

If a circulating bath is used as the water source, it may not be possible to maintain a flow rate of 200 mL per minute. While this condition will limit the lowest temperature attainable, it may be partially compensated by using colder circulating liquid. Below about -25 °C, 30% methanol in water is generally recommended as the coolant.

If the set temperature is near room temperature and too close to the temperature of the cooling water, the sample holder temperature may oscillate around the set temperature. In this case, changing the temperature of the cooling water (for example by adding ice) should stop the oscillation. Cooler circulating water can result in improved performance of the cuvette holder at low temperatures, whereas warmer water can improve performance at very high temperatures.

If a cuvette shorter than 44 mm is used, it may be difficult to remove. Use a taller cuvette instead.

4.5. Error Messages

When errors occur, the line 1 of the **TC 1** Temperature Controller display presents an error code. Line 3 of the display identifies the error and line 4 of the display presents possible solutions. The most common events that cause errors to be displayed are loose cables or inadequate coolant flow. For errors not easily solved, contact QNW at www.qnw.com/contact-us/ or service@qnw.com.

Specific error messages include the following.

5 – cell out of range

warnings: loose cable, sensor failure

The temperature controller is not receiving a reasonable response from the sensor on the cuvette tower. Either the sensor has failed or a cable is not making a good connection.

6 - cell out of range

warnings: loose cable, check cable

The temperature controller is not receiving reasonable responses from either the cell tower or heat exchanger sensors. Since it is very unlikely for both to fail, probably a cable is loose.

7 - heat exchanger error

warnings: loose cable, sensor failure

The temperature controller is not receiving a reasonable response from the sensor on the heat exchanger. Either the sensor has failed or a cable is not making a good connection.

8 - inadequate coolant

warnings: inadequate coolant, water temperature

The sensor on the heat exchanger is reading a temperature above 60 °C. Temperature control has been shut down to prevent damage to the Peltier element. Either the water was too warm or the rate of flow was inadequate to draw sufficient heat from the heat exchanger.

9 - Invalid command

The controller has been sent an invalid text command.

5. Maintenance and Spare Parts

5.1. Maintenance

The **Versa 20** and **TC 1** Temperature Controller require very little routine maintenance on the part of the user. For routine cleaning of exposed surfaces, use only a cloth dampened with water or diluted detergent. Do not use organic or abrasive solvents.

CAUTION



Caution – Any action which makes it necessary to open the **Versa 20** or **TC 1** Temperature Controller units must be executed only by QNW technicians or authorized personnel.

The water hoses and their attachments to the **Versa 20**, **TC 1** Temperature Controller, and submersible pump should be inspected prior to each usage to ensure that they are intact. Replace the nylon tubing when it becomes discolored or cracked. Periodically replace the water stored in the bucket to minimize the growth of microorganisms. When not in use, the pump, bucket, and tubing may be stored dry.

5.2. Spare Parts

Vinyl tubing and magnetic stir bars may be obtained from a variety of commercial vendors. For electrical cables to be used with **Versa 20** and **TC 1** Temperature Controller and submersible pumps, please contact QNW at service@qnw.com.

Appendix – Serial Communications for the TC 1 Temperature Controller

This document provides the serial communications protocols for version 1.0 of the firmware on the **TC 1** family of controllers.

All functions of the temperature controller can be managed from a computer, using the command set described below. If you purchased your unit as a component of a spectrometer from certain manufacturers, this feature may be implemented through traditional RS232 serial connectors on the computer or the spectrometer and on the controller. In this case they will be connected by a standard 15-pin serial extension cable (male connector on one end and female on the other). No driver installation should be needed.

Otherwise the serial linkage will be established through a USB connection between the computer and the controller. In this case the controller includes electronics which convert the USB connection to a serial communications port. However, for the port to be available to programs on the computer it will be necessary to load driver software. It is important that the driver software be loaded before connecting a USB cable between the controller and the computer. Contact Quantum Northwest at support@qnw.com for further information. Quantum Northwest can provide a control program written specifically for control of all functions of the temperature controller. Ask for program **T-App**.

In programming for the **TC 1** controller, one must adhere to the conventional notation: 8/N/1.

```

Baud:      19200
Data Bits: 8
Parity:    None
Stop Bit:  1
Flow Control:  None
  
```

For many of the commands listed below, the controller returns information in response to the command. All commands and responses are delineated by left and right square brackets ([]). Any text sent to the controller not enclosed within brackets will be ignored. In this document, an ellipsis (.....) is used to distinguish responses from commands.

Note: [command]	Purpose of the command (sent to the controller)
..... [response]	Meaning of the response (received from the controller)

1. Identify

[F1 ID ?]	What is the ID number of the sample holder being controlled?
..... [F1 ID 14]	Sample holder is a t2, Versa 20 , or other single cuvette sample holder.

The ID and version number are shown briefly on the display when the temperature controller is turned on.

Assigned Identities:

```

ID = 00 – specialty sample holder
      14 – t2
      24 – t2x2
      34 – turret6 or other multi-sample holder
  
```


- stirrer is off (+ is on, - is off)
- temperature control is on (+ is on, - is off)
- temperature is stable (S is stable, C is changing)

- [F1 IS +] Automatically report instrument status whenever it changes (e.g., due to manual changes at the controller)
- [F1 IS -] Stop automatic reports of instrument status.
- [F1 IS R] The controller has been powered off and back on again.

7. Current Sample Temperature

- [F1 CT ?] What is the current temperature of the holder?
- [F1 CT 22.84] The current temperature is 22.84 °C.
- [F1 CT +3] Periodically report current temperature every 3 seconds.
- [F1 CT -] Stop periodic current temperature reports.

8. Error Reporting

- [F1 ER ?] Report the current error.
 - [F1 ER 0] No current error.
 - [F1 ER 5] Cell T out of range (Loose cable? Sensor failure?).
 - [F1 ER 6] Cell & heat exchanger T out of range (Loose cable?).
 - [F1 ER 7] Heat exchanger T out of range (Loose cable? Sensor failure?).
 - [F1 ER 8] Inadequate coolant (check flow). Temperature control has shut down.
 - [F1 ER 9 <<bad command>>]
- Syntax error on a preceding command where <<bad command>> is the command that caused the syntax error.
- [F1 ER +] Automatically report errors when they occur.
 - [F1 ER -] Stop automatic error reports.

9. Probe Status and Temperature

- [F1 PS ?] Is there an external temperature probe connected?
- [F1 PR +] A probe is connected.
- [F1 PR -] No probe is connected.
- [F1 PS +] Enable probe status to be sent automatically when a probe is installed or removed. This is the default.
- [F1 PS -] Disable automatic sending of probe status.
- [F1 PT ?] What is the current probe temperature?
- [F1 PT +3] Periodically report the probe temperature every 3 seconds.

..... [F1 PT 22.3]	The current probe temperature is 22.3 degrees.
..... [F1 PT NA]	Probe temperature is not available.
[F1 PT -]	Stop periodic probe temperature reports.
[F1 PA S 0.5]	Set the increment for automatic reporting of the probe temperature to 0.5 degrees during a ramp. (Increment must be positive, without sign in tenths between 0.1 and 9.9 degrees, and will work for ramps going up or down.)
[F1 PA +]	Start automatic reporting of probe temperature every temperature increment (set by the command above).
..... [F1 PT 30.5]	The current probe temperature is 30.5 degrees.
[F1 PA -]	Stop automatic reporting of probe temperature every temperature increment.
[F1 PX +]	Change probe temperature returned to a precision of 0.01 degree.

10. Temperature Ramping

[F1 RR S 2.10]	Set the ramp rate to 2.10 °C/minute.
[F1 RR ?]	What is the current ramp rate?
..... [F1 RR 2.10]	Current ramp rate is 2.10 °C/minute.

For the following three commands, there are no corresponding reference commands ([R1 . . .]; see command class 13).

[F1 TL +]	Ramp the sample and reference identically.
[F1 TL -]	Ramp the sample while the reference remains stable.
[F1 TL 0]	Ramp the sample and reference independently.

The default is to ramp the sample while the reference remains stable. Ramping both the sample and reference at different rates or to different target temperatures (see below) is not possible.

To Ramp the temperature:

1. equilibrate at the starting temperature,
2. set the ramp rate,
3. set a new target temperature (command class 5).

The target temperature may be above or below the current temperature; as soon as it is set, the ramp will begin, up or down, to that new target.

After reaching the target, the controller will hold at that temperature. At any time a new ramp rate and target temperature can be set to start a new ramp.

Once you are done, set the ramp rate to 0. Otherwise, setting a new target temperature later will initiate a ramp to that target temperature.

Notes:

The minimum settable ramp rate is 0.01 °C/minute.

For higher ramp settings, the observed rate may be lower than that calculated from RT and RS or it may be nonlinear over part of the temperature range because the maximum possible rate of heating or cooling is limited (and dependent on the ramp direction as well as on the temperature).

When the ramping process is completed the controller will send a response of the form [F1 TT #] as a notification of the end of the ramp. This response can be blocked at any time by previously sending the command [F1 TT -] (see command class 5).

11. Heat Exchanger Temperature

[F1 HL ?]	What is the high temperature limit for the heat exchanger?
..... [F1 HT 60]	The heat exchanger high temperature limit is 60 °C.
[F1 HT ?]	What is the current temperature of the heat exchanger?
..... [F1 HT 39]	The current heat exchanger temperature is 39 °C.
[F1 HT +3]	Start periodic heat exchanger temperature reports every 3 seconds.
[F1 HT -]	Stop periodic heat exchanger temperature reports.

12. Cell Changing

These commands will have an effect only for a sample holder with multiple cuvette positions.

[F2 DI]	Device initialize: move to home position.
[F2 PI]	Device initialize: move to home position and reply when done.
..... [F2 DL 1]	Device is finished moving.
[F2 DL 3]	Device locate: move to position 3. (Device should be initialized prior to using this command for the first time.)
[F2 PI]	Device initialize: move to home, then to position 1 and reply when done.
..... [F2 DL 1]	Device is in position 1.
[F2 PL 4]	Device locate: move to position 4 and reply when done. (Device should be initialized prior to using this command for the first time.)
..... [F2 DL 4]	Device is now in position 4.
[F2 DD 400]	Set speed to 400 (acceptable range 100 – 900, with a default of 500).

13. Reference Cuvette

These commands will have an effect only for systems with two independently controlled sample holders.

[R1 ...]

To control and monitor the temperature and status of the reference cuvette using a Dual Temperature Controller, use any commands in classes 3-8, 10 and 11, substituting R1 for F1. There are no corresponding [R1 ...] commands for command classes 1, 2, 9 and 12.

If you wish to ramp the temperature of the reference and sample cuvettes together, please note the linking command, [F1 TL +], in command class 10.

14. Specialty Sample Holder

For a Specialty Sample Holder (ID = 00) controlled by a TC 1, there will be a second response to the [F1 ID ?] command.

[F1 ID ?] What is the ID number of the sample holder being controlled?
 [F1 ID 00] Sample holder is a **Specialty Sample Holder**
 [F1 FS (+/-)(+/-)(+/-)(+/-)(+/-)]

This response will automatically follow the [F1 ID 00] response. It defines the Functional Status (functionality) of the holder through its five (+/-) parameters. Each parameter defines whether the holder has (+) or does not have (-) the functionality as defined in the table below.

parameter #	functionality
1	Temperature Control
2	Stirring Control
3	External Probe
4	Multiple Sample Positions
5	Reference Holder

As appropriate to the functionality, the commands and responses in command classes 2 – 13 will also be available.

15. Control of Automatic Reports

The TC 1 can be controlled manually using the buttons and the display on the front. When a computer program is in use, the TC 1 will automatically send reports to the program whenever a manual change has been made. The commands in this section are intended to allow the program to prevent the TC 1 from sending those automatic reports.

[F1 XX R(+/-)] Stop (R-) or start (R+) reporting all manual changes, including those starting with [R1 and [F2.
 By replacing the XX with individual command codes, a program can selectively stop or start reporting of manual changes.

[F1 SS R-] Stop reporting manual changes for the sample stirrer.

[F1 TC R-] Stop reporting manual changes for the sample temperature control.

[F1 TT R-] Stop reporting manual changes for the sample target temperature.

[F1 PR R-] Stop reporting manual changes for the external probe connection.

[F1 RR R-] Stop reporting manual changes for the sample ramp rate.

[F2 PL R-] Stop reporting manual changes for the sample position.

[R1 SS R-] Stop reporting manual changes for the reference stirrer.

[R1 TC R-]	Stop reporting manual changes for the reference temperature control.
[R1 TT R-]	Stop reporting manual changes for the reference target temperature.
[R1 RR R-]	Stop reporting manual changes for the reference ramp rate.