

 $(\mathbf{Q}(\mathbf{Q}))$

A high-performance single-photon camera

DESCRIPTION

SPAD 512 is a photon-counting camera for high-speed imaging. The core of the camera is a SPAD image sensor with 512×512 pixels. Photon counting with up to 100'000 frames per second and zero readout noise is achieved. The global shutter enables nanosecond exposures with exposure shifts of 17 ps. The image sensor is optimized for low noise, with a typical dark count rate of less than 25 cps.

KEY BENEFITS



Widefield camera

Released in 2021, SPAD 512 is the first commercially available SPAD camera in the world. The camera features a 512×512 SPAD image sensor for both photon-counting and photon-time gating applications.



Wide detection spectra and low noise

Our single-photon detectors are fabricated in a state-of-the-art CMOS process and offer an ultra-low dark count rate of less than 25 cps. Microlenses enhance the sensor's detection efficiency.

Time gating

The sensor features time gating to study time varying signals of interest, such as FLIM. This makes it a perfect addition to any widefield FLIM microscopy setup.



Plug and Play

The system requires just a 5V power adapter and two USB3 cables to run. For full flexibility, 2 additional control lines can be connected to the SMA connectors.







APPLICATIONS

Widefield fluorescence lifetime imaging

SPAD cameras increase the overall photon throughput compared to scanned detection systems from the typical 10 Mcounts per second to 26 Gcounts per second.

Why SPAD cameras?

- Simplify FLIM setup
- Increase FLIM frame rate

High-speed imaging

SPAD cameras enable high frame rates with global shutter at zero readout noise.

Why SPAD arrays?

- Image fast phenomena in low light conditions
- Image light-in-flight





TECHNICAL SPECIFICATIONS

TYPICAL TECHNICAL SPECIFICATIONS

SENSOR	SPAD IMAGE WITH IN-PIXAL GATE
Image array	512 x 512
Pixel pitch	16.38 µm
Sensor wavelength range	400 to 900 nm
Peak photon detection probability	50% @ 520 nm
Fill factor with microlenses	>50 % for collimated light
Median dark count rate at room temperature	<25 cps
Percentage of pixels with >1 kcps	1.5%
Frame rate (max.)	100'000 fps @ 1-bit for 1 s 5'000 fps @ 4-bit semi-continuous 400 fps @ 8-bit continuous
Maximal pixel count rate with exposure modulation	167 Mcps (400 Mcps with response linearization)
Maximal pixel count rate without exposure modulation	100 kcps (240 kcps with response linearization)
Minimum gate width	6 ns
Minimum gate shift	17 ps
Exposure rise / fall time (20/80%)	350 ps / 150 ps
Lens type	C-mount



147 mm







TYPICAL PERFORMANCE CHARACTERISTICS







SYSTEM INTEGRATION

For operation, only three plugs are required, a 5 V power supply and two USB3 connections. The system software provides functionalities for photon-counting, time-gating and fluorescence lifetime imaging.

It enables 1-bit, 4-bit and 6 to 12-bit (time-gated) imaging modes and phasor FLIM processing. It can be accessed through TCP/IP for easy integration into LabVIEW, MATLAB or Python.



Pi Imaging Technology SA EPFL Innovation Park 1015 Lausanne, Switzerland info@piimaging.com www.piimaging.com +41 76 5733 314

